

Bransonic® Ultrasonic Cleaning Bath Accessories



General Description

Bransonic accessories are designed to maximize the effectiveness of your Bransonic baths. It is important that parts do not rest on the bottom of the bath, as it can inhibit the ultrasonic activity and possibly damage the unit.

A variety of perforated trays, mesh baskets, beakers, and support racks give you the option of direct or indirect cleaning. All five models come with covers to allow the cleaning bath to heat up more quickly.

Direct Cleaning Method

- Suspends parts using a perforated tray, mesh basket, or a support rack.
- Permits sound waves to penetrate directly to the parts.

Indirect Cleaning Method

- Suspends parts using a solid tray, or beakers with a beaker cover.
- Ideal for cleaning small parts.
- Discarded soil cleans easily from a beaker or solid tray.
- Indirect method is ideal for sample preparation.







Consult with your Branson representative about the Bransonic accessories and solution best suited for your application needs. **Visit us at www.Bransonic.com**

Additional Features

- Perforated trays, solid insert trays, and wire mesh baskets have built-in handles specially designed to prevent accessories from touching the tank bottom.
- Support racks allow total volume of the tank to be used while keeping parts, fixtures, or containers about ½" above the tank bottom.
- Support racks also are useful as countertop drain racks, holding parts above the countertop surface.

BRANSON

Bransonic Accessories – Order using part numbers listed under model numbers

| | Ordering Part No. | Model 1800 | Model 2800 | Model 3800 | Model 5800 | Model 8800 |
|-------------------------------------------------------------------------------------|--------------------------------------|----------------------------------------------------------------------------------------------------------------|------------------------------------------------------------|------------------------------------------------------------|---------------------------|--------------------------------|
|  | Cover (Supplied with Unit) | 100-032-515 | 100-032-517 | 100-032-519 | 100-032-510 | 000-917-050 (SS) |
|  | Perforated Tray | 100-410-160 | 100-410-162 | 100-410-164 | 100-410-166 | 100-410-168 |
|  | Mesh Basket | 100-916-333 | 100-916-334 | 100-916-335 | — | — |
|  | Solid Insert Tray | 100-410-170 | 100-410-172 | 100-410-174 | 100-410-176 | 100-410-178 |
|  | Beaker Positioning Cover | (1x600 ml) 100-032-516 | (2x600 ml) 100-032-518 (2x250 ml) 100-032-527 | (2x600 ml) 100-032-520 (3x250 ml) 100-032-528 | (4x600 ml) 100-032-511 | (6x600 ml) CPN-246-014 (SS) |
|  | Support Racks | CPN-916-039 | CPN-916-040 | CPN-916-041 | CPN-916-042 | CPN-916-043 |
|  | Beakers (All Models) | 250 ml Glass – 000-140-001 400 ml PP – 000-265-061 600 ml Glass – 000-140-004 600 ml SS – 000-410-055 | | | | |

Americas

Branson Ultrasonics Corp.
41 Eagle Road
Danbury, CT 06810, USA
T: 203-796-0400
F: 203-796-2250
Toll Free: 800-732-9262
www.bransonultrasonics.com

Europe

Branson Ultraschall
Niederlassung der Emerson
Technologies GmbH & Co. OHG
Waldstrasse 53-55
63128 Dietzenbach, Germany
T: +49-6074-497-0
F: +49-6074-497-199
www.branson.eu

Asia

Branson Ultrasonics (Shanghai) Co., Ltd.
758 Rong Le Dong Road
Song Jiang, Shanghai, PRC, 201613
T: 86-21-3781-0588
F: 86-21-5774-5100
www.branson.com.cn

All specifications subject to change without notice. All dimensions are nominal.



© Branson Ultrasonics Corporation 2012.
The Emerson logo is a trademark and service mark
of Emerson Electric Co.
Printed in the U.S.A.

BCL-0004-12